

## U-D Precision

The U-D Precision is designed for taking low volume unit dose samples from different heights within the product. The sampler takes three samples at each of three sampling heights simultaneously.


Up to 3 samples can be taken simultaneously

**Type:** Multilevel  
For Unit Dose Sampling (very low volume sizes)


**Sites:** Pharmaceutical

**Material:** 316 stainless steel

- Operation:**
1. Select the Sample Inserts of the correct volume.
  2. Insert the Sampler Inserts into the Inner Rod
  3. Push the Inner Rod into the Outer Tube so that the Sampling Inserts can not be seen through the sampling ports
  4. Insert the sampler into the product
  5. At the required depth rotate the inner core through 180° to expose the Sampling Inserts. The sample will flow into the inserts.
  6. Rotate the inner core through 180° to hide the sample cells.

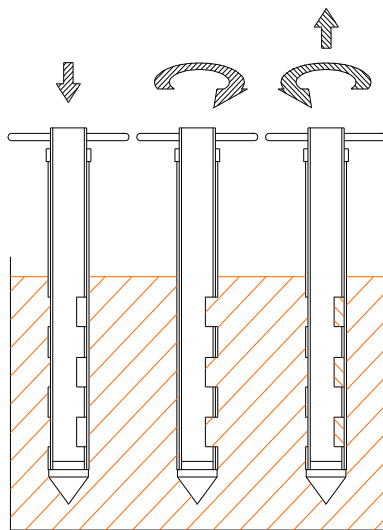

- (A) Outer Tube
- (B) Sampling Port
- (C) Inner Rod
- (D) A Set of Sampling Inserts

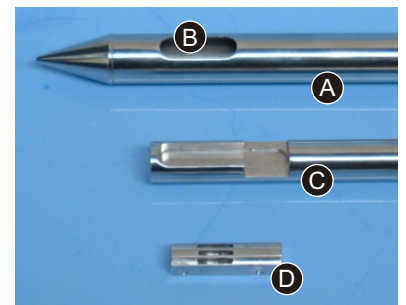

### Sampler Body

Part No.	Sampler Length	Number of Sampling Ports
I280A-600	600mm	3
I280A-1000	1000mm	3
I280A-1500	1500mm	3

### Sampling Inserts (3 inserts per Set - 3 sets will be required for each sampler)

Part No.	Volume	Price per Set		Part No.	Volume
		1 - 3 sets	4+ sets		
I280T-010	0.01ml	£148.80	£141.36	I280T-060	0.06ml
I280T-020	0.02ml	£148.80	£141.36	I280T-070	0.07ml
I280T-030	0.03ml	£148.80	£141.36	I280T-080	0.08ml
I280T-040	0.04ml	£148.80	£141.36	I280T-090	0.09ml
I280T-050	0.05ml	£148.80	£141.36	I280T-100	0.10ml
I280T-000	Blank Insert	£115.80	£110.01		